

Welcome to

BioCAM

DuplexX



User Guide Essentials

Last updated 13/08/2021

Contents

5	Introduction
6	MEA technology
6	CMOS-MEA technology
8	At a glance
9	Overview
10	HD-MEA
11	BioCAM DupleX
13	BrainWave 4 video tutorials
13	BrainWave 4
14	Precautions
15	BioCAM DupleX handling and use
16	HD-MEA handling and use
17	Cleaning procedure for HD-MEAs
18	Sterilization procedure for HD-MEAs
19	Getting started
20	Install the BioCAM DupleX
20	Setting up the BioCAM DupleX
20	Plug the HD-MEA
21	Verify proper functioning
22	Quick guide to your first use
23	Basic concepts
23	Visualization of data
24	MEA Viewer control
25	Data formats
26	Online mode (Record mode)
26	Important notes
26	Open recorder
27	Visualize live data
28	HD-MEA Calibration
29	Record live data
30	Offline mode (Playback mode)
30	Open recorded data
31	Quick analysis and save of results
32	Visualize recorded data
33	Understanding main controls affecting graphs
34	Create a group of relevant channels
35	Visualize Instantaneous MFR
36	Visualize Raster

37	Troubleshooting
38	Electrodes with no-signal (saturation)
40	HD-MEA not properly initialized
42	Not working columns
43	HD-MEA with dirty surface
44	Pseudo-Reference Electrode bubbles
45	Contact Information

1

Introduction

Welcome to the BioCAM DupleX platform and thank you for your purchase. The BioCAM DupleX platform is among the most advanced systems for managing your experiments on the emerging generation of high-density CMOS-MEAs. This manual has been written to help you in taking advantage of all functionalities provided by your BioCAM DupleX. Be sure to read this manual thoroughly, and to keep it handy when using the BioCAM DupleX platform.

Before any use of your BioCAM DupleX platform, please read the ["Precautions" P.14](#), which contains relevant information to preserve your HD-MEAs and BioCAM DupleX from possible damages.

MEA technology

Among the different methodologies used for electrophysiological measures, metal microelectrodes integrated on-chip can provide multisite measures of extracellular signals with a high signal-to-noise ratio. In addition, by applying voltage or current stimuli to the same microelectrodes it is possible to depolarize cells or tissues, thus establishing bi-directional interfaces. As established over several decades of research, both sensing and actuating performances of microelectrodes can be applied to study a wide range of electrogenic cells and tissues, including neuronal and cardiac preparations.

Conventional microelectrode arrays (MEAs) are bio-sensing chips realized by means of thin-film technologies and do not integrate any on-chip microelectronic circuit. Therefore, conventional MEAs are passive devices made on silicon, glass or polymeric substrates. Each microelectrode can be made by different materials (e.g. Pt, IrOx, TiN) and it is individually wired on-chip and connected to an external amplifier and data acquisition (DAQ) instrument. Due to the need of this MEA technology of individually routing on-chip each electrode, space constraints and wiring encumbrance impede the realization of dense and large electrode arrays. Thus, for conventional MEAs the typical electrode pitch is in the range of 100 μm and the array includes from 60 up to 256 microelectrodes. In addition, given the distance between the electrodes and the off-chip amplifiers and the use of interconnecting wires, conventional MEAs are subjected to inductive coupling noise.

CMOS-MEA technology

The electrode density and array sizes can be increased by changing the technology used to realize microelectrode arrays (MEAs). High-density MEAs are realized with complementary metal-oxide-semiconductor technology (CMOS), as it is done for microelectronic devices (e.g., computer microprocessors) and light-imaging devices (e.g., camera sensor), and with post-processing methods to optimize the electrode performances.

In short, the CMOS technology allows the realization of active electrode-pixels integrated in small areas of a few square micrometers electrodes, amplifiers and signal conditioning circuits in-pixel, just underneath each electrode site. On-chip, additional amplification stages, time-division multiplexing and high-speed addressing circuits are provided.

In particular, the circuit architecture of 3Brain's HD-MEAs is based on the Active Pixel Sensor concepts, as commonly used for light imaging CMOS cameras, and was therefore designed to allow full array recordings at sufficiently high sampling frequency for each electrode.

In HD-MEA, the local in-pixel buffers underneath each electrode allow to locally adapt signals from high impedance (electrode-side) to low impedance (wiring-side) to avoid the induction of coupling noise from electrical wiring before amplification. Additionally, the on-chip multiplexing and addressing circuits allow the minimization of the number of wiring outputs despite the large number of electrodes that is to be simultaneously measured. Indeed, each electrode signal is not individually wired off-chip. Rather, signals of multiple electrodes are multiplexed at high frequency over a few number of wires. In this way, several thousands of electrodes can be recorded on only a few tens of output wires.

The in-pixel circuit implemented in HD-MEAs allows recordings from a large signal bandwidth that includes field potentials and spikes. Instead of implementing an AC-coupling solution that would require large areas for the integration of adequate capacitors, the HD-MEAs in-pixel circuit integrates an auto-zeroing circuit that is regularly calibrated to the electrode DC voltage and that subtracts this calibrated DC voltage from the electrode signal before the first amplification stage. This process is called calibration and it also allows the adaptation of the circuit performance to different experimental conditions. In particular, all CMOS devices are photosensitive and photo-generated charges give rise to different DC drifting times that can bring in-pixel amplifiers to saturate. HD-MEAs mitigate these effects and hence allow experiments under light stimulation conditions (as required for instance for retina stimulation) using higher calibration frequencies to keep the DC input signal at the working point of the amplifier and to avoid the saturation of the amplifier.



2

At a glance

Overview

The BioCAM DupleX is the latest generation of our single HD-MEA recording platforms. Similarly to the BioCAM X, it allows the acquisition of data from our CMOS high-density MEAs. However, with BioCAM DupleX's improved and completely re-designed electronics, this provides an advanced on-line data-processing compatibility with the latest Accura MEAs. The Accura MEAs feature recording/stimulation from any of the 4096 electrodes, an integrated circuitry to handle the reference ground on chip and a motorized locking mechanism to improve connection reliability of the system with the MEA.

Either your biological models are primarily neuronal cultures, human-derived stem cell preparations, cardiac cells or brain tissues as slices, retinas or organoids, the optimized electronics of the BioCAM DupleX will be able to detect both slow oscillations (field potentials) as well as spiking activity with unprecedented sensitivity. Examples of recordings provided by different laboratories using the BioCAM platforms are available on the 3Brain website and uploads of novel examples based on your research are warmly welcomed.



Overall, the platform consists of three basic components:

1. A multiuse HD-MEA cartridge available in different models and incorporating the CMOS-MEA silicon chip;
2. The acquisition BioCAM DupleX hardware that reads out electrophysiological signals from the HD-MEA; and provide electrical stimulation through the electrodes (Accura chip model only).
3. A high performance host PC with pre-installed BrainWave 4, our latest all-in-one software generation to acquire, visualize, store and analyse the data recorded by the BioCAM DupleX.

HD-MEA

The HD-MEA cartridge integrates on an electronic substrate, the CMOS-MEA chip and a reservoir chamber to maintain cultures or tissues that goes in contact with the electrode array either under cell culture media or perfused media, respectively. The cartridge is 54 mm x 54 mm (2-1/8 in x 2-1/8 in) in size and is conceived to be placed in an incubator for cells or tissue cultures.

- *Contact pads* : gold-plated pads for connection between the HD-MEA and the BioCAM Duplex. Properly cleaned and non-oxidized pads are needed for a stable connection.
- *PCB* : the printed circuit board substrate of the HD-MEA. Different colors are available and according to the layout of the CMOS-MEA.
- *Reservoir* : a bio-compatible plastic chamber for the bath solution used for the biological preparation. The volume of solution that can be contained is of about 2mL. The external ring defining the reservoir volume has an external diameter 28mm (1-7/64 in), internal diameter 25mm (0-63/64 in), height 5mm (0-13/64 in).
- *CMOS-MEA* : chip integrating thousands of active electrodes for high-resolution electrophysiological recordings from cells or tissues placed on its active area. Different application-specific MEA layouts according to the HD-MEA model are available.
- *Reference Electrode* : two large gold pads integrated into the reservoir to provide the reference ground. Implemented to avoid the usage of an external ground reference.

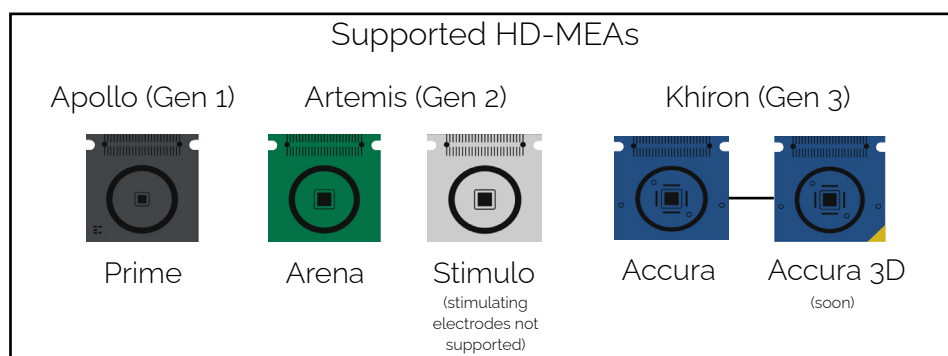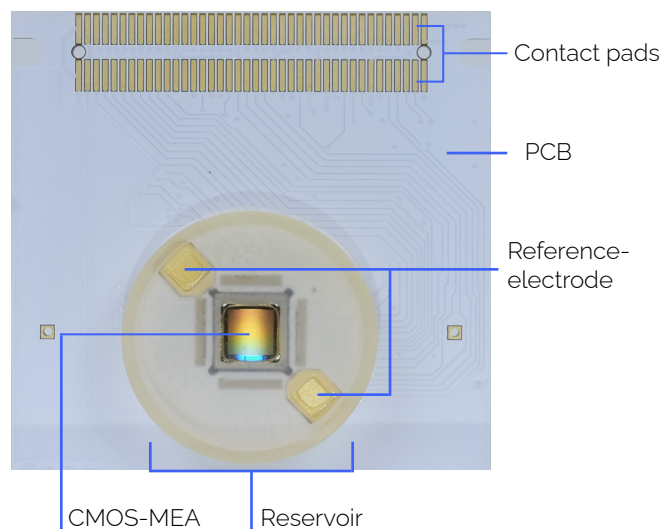


BioCAM Duplex


The BioCAM Duplex is the station where HD-MEAs are plugged into for electrophysiological measures. This instrument is compatible with all MEA Probe types including also those with electrical stimulation capabilities. It contains all the electronics, real-time hardware and logic to control the CMOS-MEAs chips on the HD-MEA cartridges and to acquire and pre-process large-volume of electrophysiological data before sending it to the host PC.



- *Lock button* : automatically locks the HD-MEA chip into the bay when initially pressed. Pressing again will release the chip. Lock/unlock status can be checked by the LED lock light indications.
- *LED light display* : BioCAM Duplex provides a set of LEDs to monitor powering, recording and chip lock/unlock status
- *Peltier-Element* : active element to control the temperature of your preparation on HD-MEAs
- *HD-MEA bay* : bay for inserting the HD-MEA and connecting it to the BioCAM Duplex.

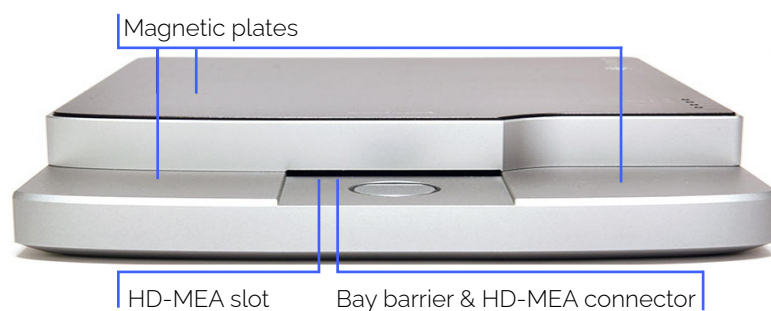

- *Magnetic plates* : ferromagnetic stainless steel to attach magnetic holders (e.g., for perfusion).
- *HD-MEA slot* : the slot through which the HD-MEA (contact pads side) is inserted into the Duplex for electrical interfacing.
- *Bay Barrier & HD-MEA connector* : immediately behind the HD-MEA slot, a protective movable Bay Barrier blocks accidental overflows of liquids that might get in touch and damage the internal electronics. Behind the Bay Barrier, the HD-MEA connector allows to connect the HD-MEAs contact pads to the internal circuitry. Electrical contacts are established only once the locking system is activated.

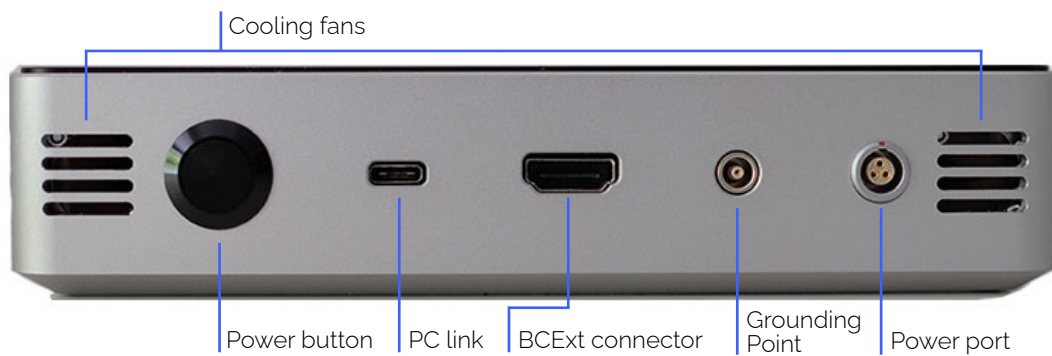

- *Cooling fans* : releases excess heat generated from the core of the Duplex. It is mandatory to allow airflow and not be blocked directly against any obstacles.
- *Power button* : push-button to power ON or OFF off the Duplex (the instrument needs to be connected to a power line using the provided adaptor).
- *PC link* : the UCB-C provides a connection for streaming the electrophysiological data from the Duplex to the host PC. It is also used by BrainWave 4 to provide BioCAM Duplex firmware updates.
- *BCExt connector* : plug the external connector box (BX110A) to access extended functionalities such as triggering and stimulation.
- *Power port* : plug for the power supply unit.


BrainWave 4 Video Tutorials

Scan this QR Code to get to multiple BrainWave 4 Video tutorials that we've set up for additional guidance. Direct link is : https://www.youtube.com/user/3BrainGmbH/playlists?view=50&sort=dd&shelf_id=1



BrainWave 4

BrainWave 4 is a complete software application developed to exploit all the functionalities of your Bio-CAM Duplex platform. It is equipped with advanced tools adapted to manage high-resolution electrophysiological data acquired from HD-MEAs. When launched it appears as follows (without experiments ongoing):

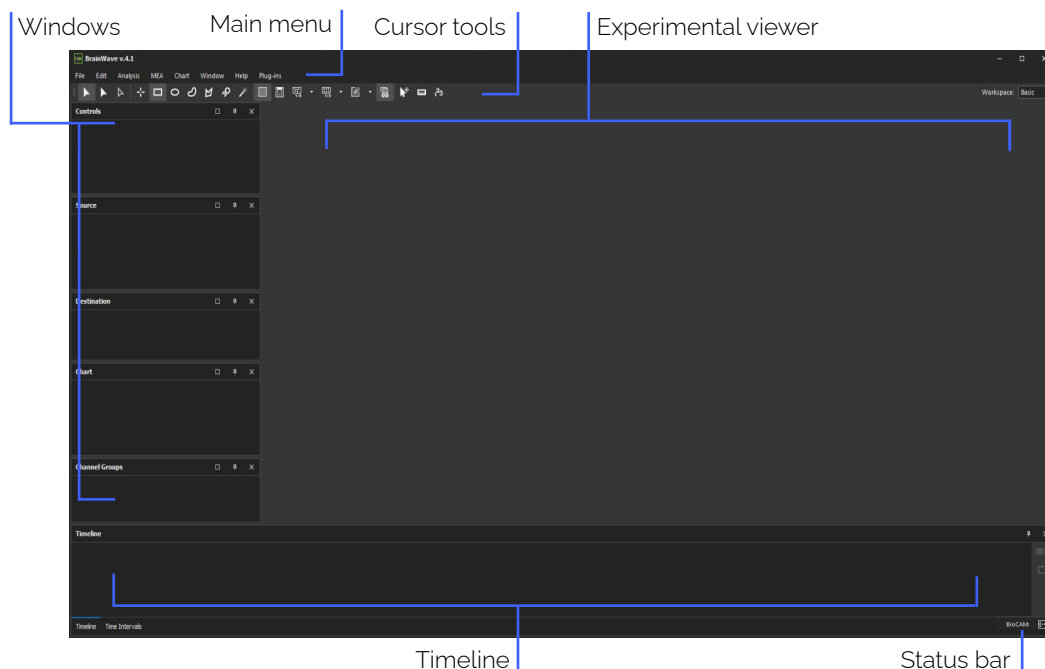

- *Main Menu*: access main functions by clicking on the main menu.
- *Cursor tools*: here you can change the behavior of your cursor
- *Windows*: contains dockable windows that are automatically updated according to the currently selected experiment loaded into the Experimental viewer. Windows allow you to modify the experiment's options and parameters which can allow you to modify different properties of your experiments. These windows can be shown or hidden by choosing the Window menu on the Main Menu
- *Experimental Viewer*: arena for visualizing online and offline experiments through tabs.
- *Right Status Bar*: notifies you about the status of the controlled devices. It displays LED icons for Bio-CAM Duplex and HD-MEA devices. If the LED icon on the left side displays with "DX" in representation of the Duplex, then it is properly connected and recognized. The display on the right "[...]" will light up when a HD-MEA is correctly placed and locked in as well.

[...] Disconnected

DX Duplex
Recognized

DX Duplex & HD-MEA
Recognized

DX ! Duplex or HD-MEA
Error (hover for details)



3

Precautions

BioCAM DupleX handling and use

The BioCAM DupleX instrument is made with a solid metal case and with optimized electronics that make the system robust to mechanical and vibrational noise. Even if it is recommended, the BioCAM DupleX's installation on an anti-vibration table is not mandatory.

Moreover, the instrument has been design to be shielded from external electrical environmental noise. However, to guarantee optimal recording performances simple precautions as listed below here should be followed:

- Position the BioCAM DupleX far from potential strong noise sources, such as fridges, incubators etc... Indeed, if not well shielded these instruments might generate electromagnetic noise.
- Connect the BioCAM DupleX to adequate power lines with low noise levels.
- Take care in avoiding that the electrical cables connected to the BioCAM DupleX (in particular, the one for the power supply and for the PC Link) do not pass near other power supplies.
- Check that the cable connectors are well inserted into the BioCAM DupleX and the host PC before operating the platform.

The system allows for the integration of a perfusion system as it is needed for maintaining tissues during electrophysiological measures. Even if precautions have been taken in designing the system to prevent liquids in entering into contact with the internal electronics, spill outs of electrophysiological solution that might get into the HD-MEA slot should be strictly avoided. However, should this happen please follow the following steps:

1. **Immediately unlock and remove the HD-MEA and switch OFF the BioCAM DupleX.**
2. Accurately clean the HD-MEA Bay of the BioCAM DupleX with ethanol >96% and let it dry. Execute the same procedure for all the HD-MEA parts but for the HD-MEA reservoir and in particular for the HD-MEA connector pads.
3. **In case after the cleaning with ethanol, there are still doubts that some liquid could have penetrated into the BioCAM DupleX, do not switch on the system for at least one day, allowing the liquid to be completely dried and preserving internal circuitry from potential oxidative and short-circuit effects.**

HD-MEA handling and use

HD-MEA are low-power active electronic circuits and it is highly recommended to avoid the following operations that might result in a damaging of their electronics:

- Do not touch the contact pads of the HD-MEA. Electrostatic charges might damage the on-chip circuits and might result in damaging the device. Always handle the HD-MEA by holding it on the sides and ideally by using plastic gloves.
- Do not hold the HD-MEA by its reservoir to avoid detaching it.
- Do not touch the active area of the HD-MEAs with any tools except those specifically allowed (see ["Cleaning procedure for HD-MEA" P17](#)). In particular, hard or semi-hard objects (e.g. metal or plastic tools) can irreparably damage the electrodes and the on-chip circuits.
- For cleaning or for sterilization, avoid immersing the entire HD-MEA in water or ethanol. Prolonged immersion in water of the entire HD-MEA might cause oxidation of the contact pads. In ethanol, some parts of the HD-MEA might be deteriorated. As a general rule only the chamber can be wet while the rest of the chip should stay dry.
- It is highly suggested not to place the HD-MEAs in autoclaves, ovens or under UV-light exposure for sterilization. These methods might deteriorate the CMOS devices, resulting in malfunctioning or in reduced lifetime. For the HD-MEA sterilization, please refer to the ["Sterilization procedure for HD-MEAs" P18](#).
- Maintain the pH of the electrophysiological solutions used for neuronal cultures or brain tissues possibly at physiological conditions (7-7.5). Important changes in the pH of the solution might damage the electrodes.

During recordings (both cultures and brain tissues) users have to comply with the following precautions:

- Before contacting the HD-MEA into the BioCAM Duplex, clean the metal pad contacts with a tissue paper soaked with ethanol 96% and let it dry for few seconds.
- Avoid liquids to spill out and strictly avoid liquids in entering into contact with the pads (the BioCAM system is equipped with a Bay Barrier to protect the connecting socket). In case users experience liquid spill out, follows instructions provided in the ["BioCAM Duplex handling and use" P16](#) section.

In order to ensure a good quality of the recordings:

- Avoid using HEPES (4-(2-hydroxyethyl)-1-piperazineethanesulfonic acid) in the solutions used for electrophysiological recordings, since we observed that it can interfere with the electrodes, making the chip unstable and the recordings very noisy.
- HD-MEAs are light-sensitive devices and their performances might be affected by a too highly intensified direct exposure or by noisy light sources.

Cleaning procedure for HD-MEAs

Normal cleaning

It is a good practice to clean the HD-MEAs immediately after the experimental recordings both for cell cultures as well as brain tissues. First, rinse the HD-MEAs abundantly with Double Deionized Water (DDW), then fill the chambers with a detergent such as WPI-Enzol (WPI) or Terg-A-zyme (Alconox) and gently pipette for a few seconds. Leave the detergent for a few minutes (typically 3-5 min) and then pipette again.

After the use of the detergent, rinse the HD-MEAs abundantly with DDW, then leave the chamber filled with DDW for 1-2 minutes and repeat this operation 3-4 times in order to be assured of washing out the detergent completely.

Intense cleaning

In case the chips are very dirty, users can mechanically remove cellular debris with a soft brush. Be aware that this operation can damage the sensing area of the chip, so it is suggested only if, after a normal cleaning procedure, the chips still presents dirtiness and in any case it has to be performed very mildly and carefully.

Finishing and storage

HD-MEAs can be dried by using a gentle flux of Nitrogen air (do not directly expose the recording area to the flux to avoid potential electrode damages). Alternatively let the HD-MEAs dry on a bench by covering them with a plastic petri dish to avoid dust deposition on the recording area.

The HD-MEA's area out of the chamber can be cleaned with a tissue soaked in ethanol >96%. It's a good practice to always wear gloves while handling the chips.

Once the HD-MEAs are dried they should be stored in a closed box in order to protect them from dust and dirtiness.

As a general rule, avoid sterilizing or using ethanol in the HD-MEA chamber without having preventively cleaned the chip from cellular debris of previous experiments. Alcohol can fix biological materials on the recording area of the chips, however it will result in a degeneration of the signal recording performances.

Sterilization procedure for HD-MEAs

As indicated in the previous section, it is highly suggested not to place HD-MEAs in autoclaves, ovens or under UV-light exposure for sterilization purposes. These procedures risk deteriorating the chips resulting in not optimal recording performances and a shorter chip lifespan.

To sterilize chips, first perform a cleaning procedure as indicated in the previous ["Cleaning procedure for HD-MEAs" P.17](#) and adopt the following procedures (that has to be performed under a biological hood):

1. Clean the areas of the HD-MEA outside the reservoir (including the external part of the reservoir itself) by using a tissue soaked with ethanol 96% and put it inside sterile plastic petri dishes.
2. Fill the reservoir completely with ethanol 70%. Also dip the border of the reservoir (the tiny upward-oriented side of the ring) with the same solution. Proceed to wait for 20 minutes.
3. Suck the ethanol with a pipette or with a vacuum pump, always using sterile tips.
4. Fill 3/4 of the reservoir of the HD-MEAs with sterile DDW, wait a few seconds and always suck the liquid with sterile tips. Repeat these operations 4 times in order to ensure of washing out the ethanol completely. When removing sterile DDW for the last time, be sure to have completely dried the chamber. Aspirate the liquid on the recording area by always keeping the tip close to the reservoir/CMOS border but not touching the CMOS to avoid any risk in damaging the electrodes.



4

Getting started

Installing BrainWave 4

The Brainwave 4 software will be found already installed with your BioCAM Duplex platform. However, due to any other circumstances you can still search, download and install BrainWave 4 directly from this link: <http://www.3brain.com/resources/downloads>

Setting up the BioCAM Duplex

1. Plug the power supply cable into the power port located in the back of BioCAM Duplex and then connect the power supply into an electrical outlet. Switch ON the Duplex by pressing the power button on the back. The Duplex's Power LED light should turn ON.
2. Turn on the host PC, plug one end of the PC link into the USB C port on your host PC and the other end of the PC link to the Duplex. The USB Type-C (v.3.1) cable is provided with the Duplex. If you require another, please ensure to purchase the Type C with at least or above version 3.1.
3. Launch BrainWave 4 and verify that the Duplex is properly recognized by checking that the LED icon in the BrainWave's Right Status Bar is turned ON, with the display of DX to indicate it's recognition of the Duplex platform, as indicated in previous section ["BrainWave 4" P.13](#).

Plug the HD-MEA

HD-MEAs are connected to the BioCAM Duplex through the lock system. Once a HD-MEA is inserted into the HD-MEA bay of the BioCAM Duplex, press the lock button to initiate the locking mechanism.

1. Insert the HD-MEA into the Duplex sliding it along the HD-MEA bay. The HD-MEA should enter into the Duplex's HD-MEA slot till the distance between the end of the HD-MEA and the Duplex's end is approximately about 8mm(1/3 in). If the HD-MEA is not at it's end position, the LED for the locking indication will display red. Repress the lock button to release the lock and try to re-insert the HD-MEA again




2. Once you've push the lock button, and the locking mechanism has recognized the correct placement of the inserted HD-MEA. The LED lightning indicators on the device should turn ON, as indicated on the next section ["Verifying proper functioning" P.21](#) . BrainWave 4 should recognize the input of the HD-MEA with it's Status Bar displays, as shown on ["BrainWave 4" P.13](#).

Verify proper functioning

Once you have installed and turned on the BioCAM Duplex, in case there are any updates, BrainWave will notify you to update on the host PC. You can easily verify the proper functioning of the system by the following check list:

1. Verify that both Duplex LED and HD-MEA LED icons in the BrainWave's Right Status Bar are ON (see section ["BrainWave 4" P.13](#)).
2. On BrainWave go to File > New Recording Session. On the new opening window click on the Play button to start, the "Status" LED on the BioCAM Duplex will turn on. Select a few channels to open their graphs (see ["Visualize live data" P.27](#)) and verify that the data is acquired from the HD-MEA (if you have not placed a living sample on it, you should visualize noise traces in the range of +/- 40 μ Volt)

If anything of the above went wrong, please check ["Troubleshoot" P.37](#). If the problem is still not solved, contact the customer support.

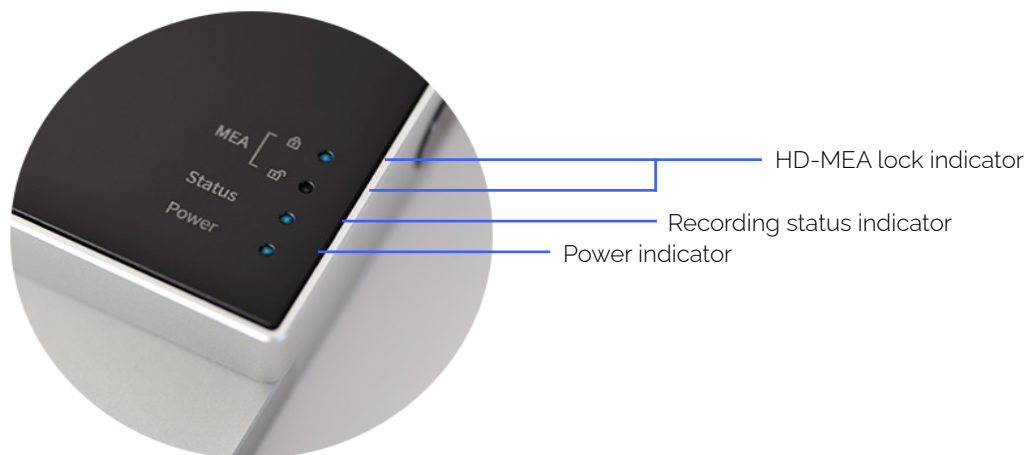

Unplug the HD-MEA :

1. Press the BioCAM Duplex's lock button, blue lights should move from the locked to unlocked on the HD-MEA lock LED indicator.
2. Once the light indicates the HD-MEA as unlocked, simply remove from the slot.



5

Quick guide to
your first use


Basic concepts

Visualization of data

BrainWave 4 makes use of tools designed to manage high resolution data coming from CMOS-MEAs. This has been implemented by combining conventional approaches with innovative ones. In particular, since HD-MEAs act as "cameras for extracellular electrophysiological potentials", several functionalities were introduced following similar approaches as used in multimedia.

BrainWave 4 is an all-in-one software including tools for recording, visualizing, analyzing and exporting your data. Use of BrainWave 4 differs based on whether you are visualizing live data from your biological preparation or whether you are playing back data previously recorded. These two different scenarios are referred as online mode and offline mode respectively.

The term Experiment can be used both for an online and an offline experiment. When a new Experiment is started (online) or opened (offline), its data is loaded and displayed in the BrainWave 4's Experimental viewer area. More Experiments can be loaded during a single application session, in particular you can manage at the same time multiple offline Experiments and one single online Experiment. When multiple Experiments are open, they are organized in tabbed documents. When passing from an Experiment to another by clicking on its corresponding tab, BrainWave 4 updates all the windows with the Experiment's data.

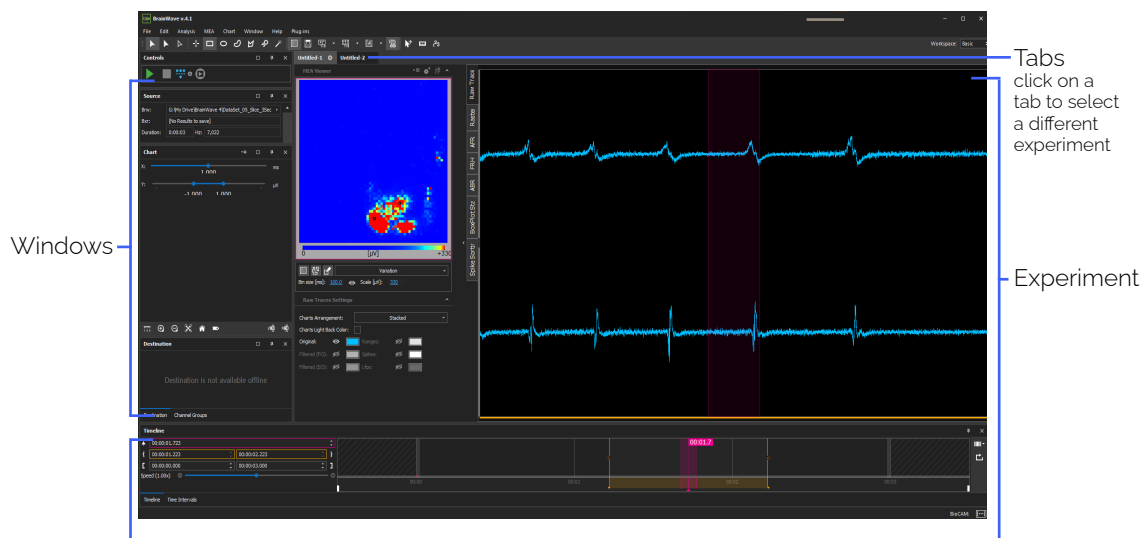

In Brainwave 4 there are two experimental visualization modes.


Online : Real time, when you access the live data from a biological sample on a HD-MEA chip. This data can be visualized directly and/or recorded to your hard disk.

Offline: refers to playback and analysis of recorded experiment's data

MEA Viewer control

Electrodes integrated in the CMOS-MEA are referred in BrainWave as Channels (abbreviated Ch). A fundamental control to browse your data is the MEA Viewer

MEA Viewer displays the activity at the current time position by a false-color map. The map gives a representation of the activity by converting the channels' voltage values in your MEA to pixel colors. This is done according to the color scale below and settings. At a specific time location in your recorded Experiment the MEA Viewer gives an image representation of the activity. When instead you visualize live data acquired from the HD-MEA or when you playback recorded Experiments, the MEA Viewer refreshes continuously, thus creating an activity movie.

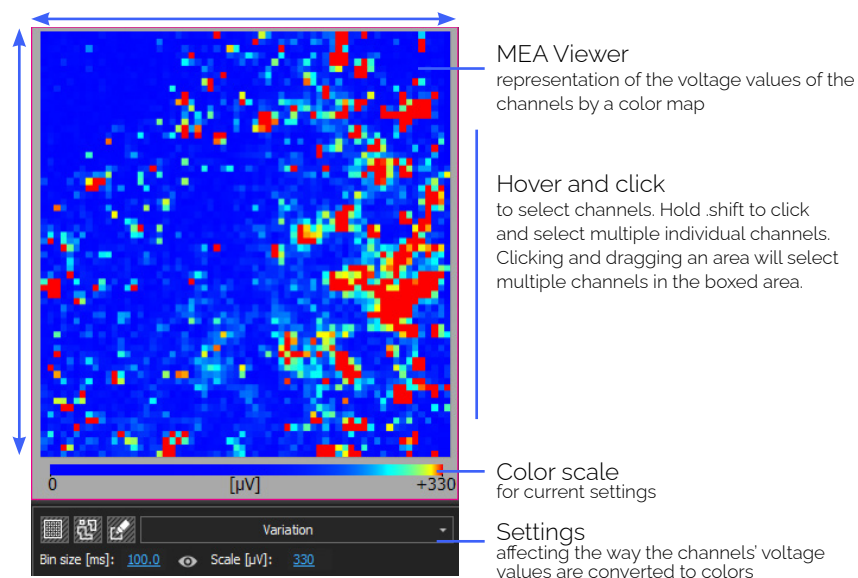

Each square pixel in the MEA Viewer represents a channel in your MEA. Channels are identified by their row and column coordinates. By convention row coordinate comes first than column coordinate, hence for instance the channel [10, 5] is the channel at row 10 and column 5.

MEA Viewer allows you to select single or multiple channels. To select multiple channels you can either hold down Shift on your keyboard while clicking on channels or click and drag an area over the MEA Viewer.

Data formats

BrainWave 4 works with two main file types although it can export data in other formats. Those two own file types are BRW- and BXR-files. Both file types are using HDF5 hierarchical data format (hdfgroup.org). You can easily open them within most operating systems (including Windows, Linux and Mac OS) and most used data analysis environments, such as, among others, Python, Matlab, Scilab, Octave and R. You can find detailed documentation on both data files on 3brain.com/resources/downloads.



BRW stands for Brainwave RaW data file and contains continuous or quasi-continuous data streamed from the the BioCAM. BRW files, according to your acquisition settings, can contain for instance raw data recorded from all the HD-MEA channels or a temporal subset of those data. The following section will teach you how to record plain raw data file.



BXR stands for Brainwave eXperiment Result data file and contains discrete data obtained by analyzing and processing the raw data. BXR files can for instance, contain the timestamps of the detected spikes. In the following section you will learn to create a BXR-file starting from a BRW-file.

Conceptually, a BXR-file is created starting from a BRW-file, hence a BXR-file is associated and linked to its parent BRW-file (while it is not true that a BRW-file is linked to a BXR-file). BXR-file can be deleted and recreated starting from the BRW-file, while if a BRW-file is deleted the contained data is lost forever. However, since BRW-file are typically large files containing a huge amount of information, sometimes you might find useful to delete the BRW-file once the relevant results have been obtained and saved in a BXR-file.

Online mode (Record mode)

Important notes


Before moving forward, follow all the steps detailed in ["Install the BioCAM Duplex" P.20](#) and ["Plug the HD-MEA" P.20](#). Hence, make sure that the BioCAM Duplex and the HD-MEA are properly connected by checking that the corresponding LED icons in the BrainWave 4's Right Status Bar are turned ON (blue icon).

Open recorder

Open the recorder by either

- Choosing File > New Recording Session, or
- Right-clicking on the Workspace > New Recording Session, or
- Pressing Ctrl+N

A new online Experiment opens as a tabbed document. Only one online Experiment can be open during an application session and it can be identified by the icon and the "Recorder" text on the document tab.

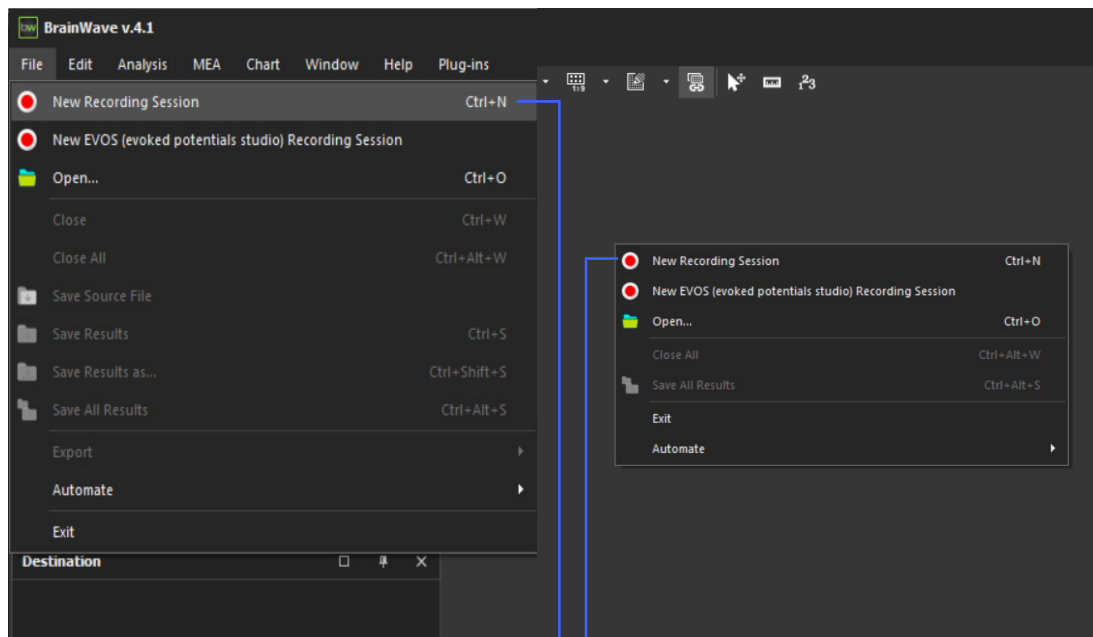

File > New Recording Session
or Pressing Ctrl+N

Right-clicking on the Workspace >
New Recording Session

Visualize live data


To start an acquisition and visualizing data read out from the HD-MEA:

1. Click Play under the Commands window to start the acquisition. The MEA Viewer starts refreshing and displaying data.
2. Click on the MEA Viewer to select individual channels or Shift-click to select more channels. For each selected channel a new chart is plotted on the right showing the voltage data being read out from that channel.
3. Use the Chart window to adjust channels' charts. Two controls are available on the Chart window for both the time (X) and amplitude (Y) axis scale values of the charts. The time control regulates the time window used by the charts. By selecting a given time window width (e.g., 500ms) the charts are refreshed by periods of the same time duration. The amplitude control allows the selection of the Minimum and Maximum values visible on the charts in the Y axis.

To stop the acquisition simply click on the Stop button under the Commands window.

1. Click Play to start an acquisition.

MEA Viewer starts to display data coming from the HD-MEA by converting MEA channels' voltage values into a colored movie.

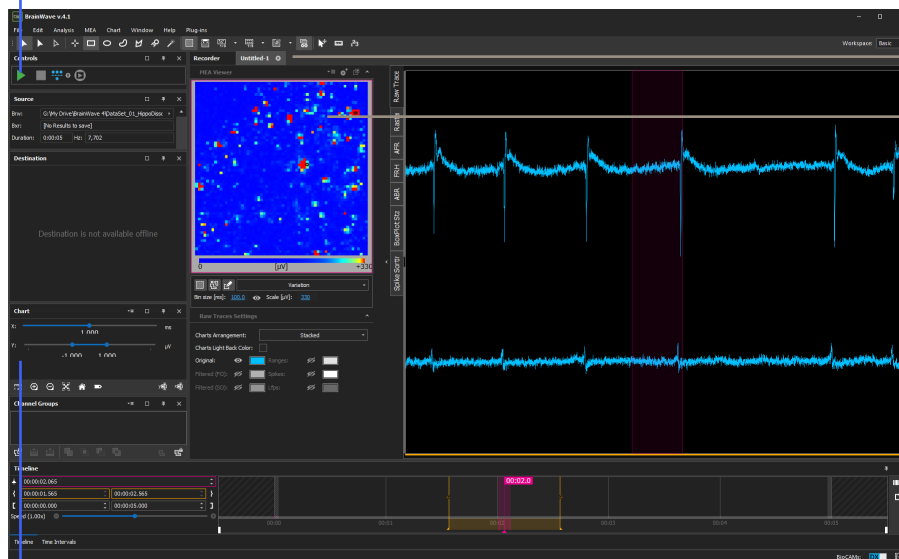

Recording tab

2. Select Channels by clicking on the MEA Viewer. To select more channels hold Shift.


Selected channels are visualized in individual charts.

3. Chart window to adjust charts scale by dragging bars for time window width and for amplitude range.

HD-MEA Calibration


In the case of when you experience several electrodes within short or prolonged periods of time with no visible signals, you may be experiencing a signal saturation problem. In this case, follow these steps before proceeding on with recording data.

A typical example of a HD-MEA with saturating behavior is illustrated in the image below. On the left, the MEA Viewer shows many deep blue colored pixels indicating not working channels while on the right typical raw data traces with short periods of saturation (top-left, top-right and bottom-left) or even with persistent saturation (bottom-right) are shown.

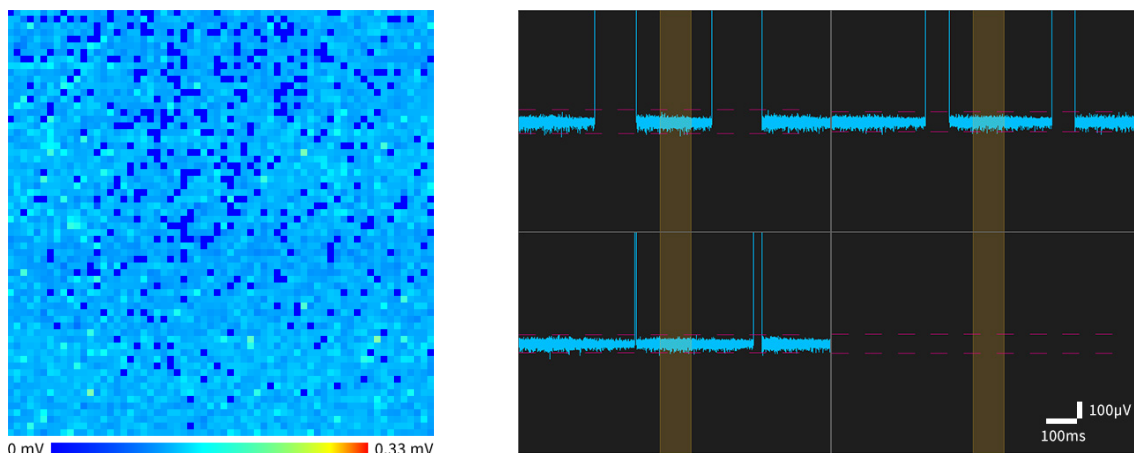

This behavior might be related to a wrong setting of the working point of the amplifiers that are integrated underneath the electrodes. In such case, follow the instructions below for a correct HD-MEA calibration; if the problem persists you can refer to the section ["Electrodes with no-signal \(saturation\)" P.38](#).

BrainWave features two functions to perform HD-MEA calibration: an automatic and a manual procedure. The automatic calibration button is positioned on the right of the window. To perform an automatic calibration you first need to start an acquisition (by clicking on the Play button) and then to click on the Auto button. The procedure will take from few tens of seconds up to roughly a minute, depending on the HD-MEA conditions.

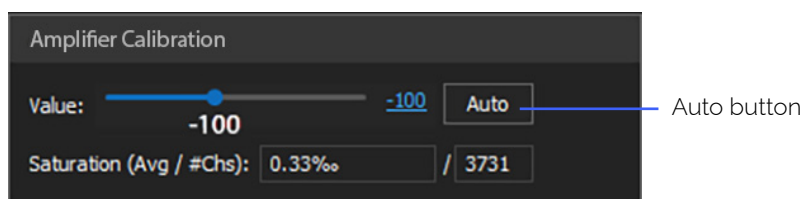


If automatic calibration is not satisfactory you can perform a manual calibration. From the window you can move the sliding bar to set a value for the amplifier bias that minimizes the saturation of the electrodes. Normally, value in the range of Prime Arena and Stimulo: -300 to -400 and Accura: 0 to 150 allows to achieve good performances, but values out of this range can also be needed. Moving the bar toward higher negative values will eliminate saturation from those electrodes that oscillate between working condition to saturation (i.e. blinking deep blue pixels in the MEA Viewer control). If you set a too high negative value, some electrodes will start saturating permanently (i.e. fixed deep blue pixels in the MEA Viewer control). For a correct setting start from values close to 0 and move the bar by small steps towards the left until an optimal condition is found.

Record live data

Complete the steps to ["Visualize live data" P27](#) and then:

4. Click the browse button on the Info window to choose the path where to save your data. Be aware that to safely record from high resolution MEA you need high performance hard disks. Hence, browse to a file located either on a high-end disk (e.g., SSD; [solid-state drive](#)) or on a disk pool configured with RAID 0 ([striping](#)). If your order included a pre-configured host PC, the disk partition where you should record your file is typically labeled with LiveRec or RAID0. Once you chose the path, by default Brain-Wave 4 selects for you the file name, but you can modify this by typing in the Name field.
5. Choose the length of the experiment you want to record by bearing in mind that the experiment will last this duration or shorter, in case you stop the recording ahead of time, but not longer.
6. Click the Play button in the Commands window to start recording. In the Progress window you can see the elapsed and remaining time of the recording.
7. Wait until the set experiment length or click the Stop button at any time to finish recording. The recorded file is a BRW-file located at the position you selected in step 4.

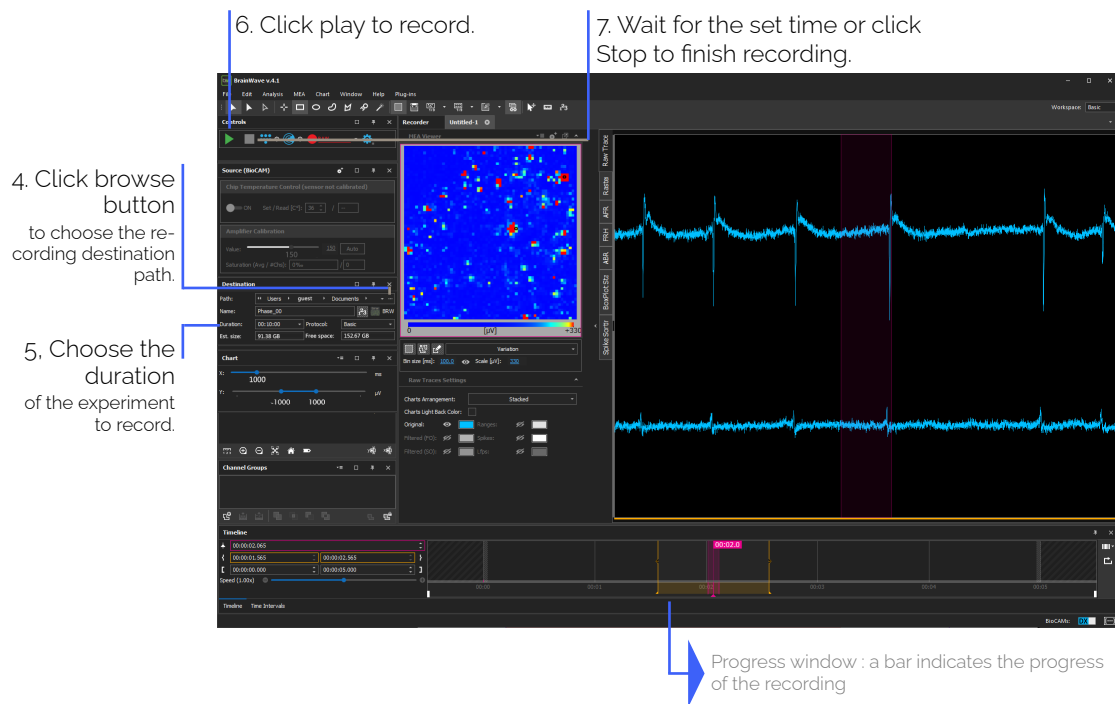

Offline mode (Playback mode)

The tutorial contained in this section is based on the demo file on dissociated neuronal cultures you can find on 3brain.com/resources/downloads. Download that file and then extract it to a known location by using 7-Zip or WinRAR.


Open recorded data

Locate the demo BRW-file to open. Then to open it in BrainWave 4 do one of the following:

- Choose File > Open, or
- right-click on the Workspace > Open, or
- press Ctrl+O, or
- drag the file from Windows Explorer into BrainWave 4, or
- double-click on the file in Windows Explorer.

A new offline Experiment opens as a tabbed document. The tab name is the name of the BXR-file associated to the BRW-file. Since you have open a BRW-file and no BXR-file has still been created the tab name will be in the form "Untitled-X", being X a progressive application session-wide number. Hence, if this is the first file you open in this session it will be named "Untitled-1".

The offline Experiment tab and the windows are similar to the online case, however some controls and functions change.

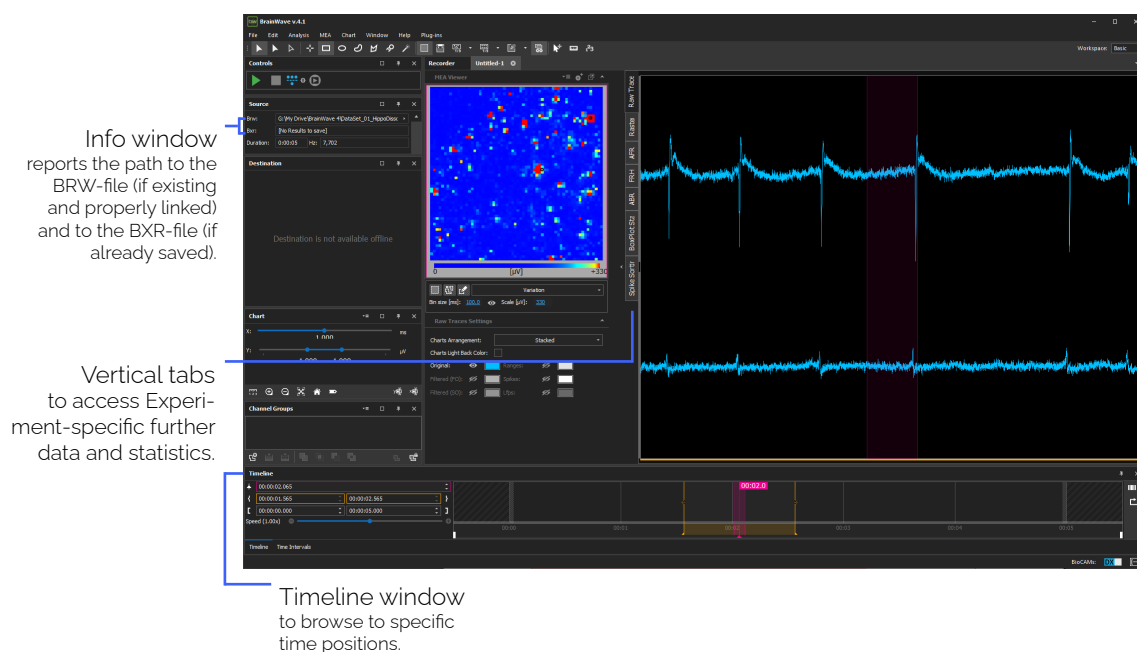

Quick analysis and save of results

As a first step, we perform a spike detection on the loaded data. Choose Analysis > Peak Detection from the Main Menu or alternatively press Ctrl+P. On the new opening window, simply leave the default parameters and click OK and then wait for the analysis to be accomplished.

Once finished, the BXR-field in the Info window turns yellow to indicate that there are unsaved data. To save data do one of the following:

- Click on the icon right to the BXR field in the Info window;
- Click File > Save
- Press Ctrl+S


Once you have saved the data, the Experiment tab name indicates that the BXR-file name and the BXR field in the Info window are updated too.

Visualize recorded data

As a consequence of the analysis, notice that the Timeline bar is now filled with vertical lines whose colors indicate the degree of the overall detected activity (averaged on the entire MEA) along the temporal axis. A blue-color indicates a relatively low activity, while a red-color a relatively high activity around that time position.


To browse your Experiment for signals:

1. Click the MEA Viewer on channel [24, g] to open its chart
2. Click the Timeline bar in correspondence of the first red vertical line (near 0.2s). You will see the activity around the selected time position for the entire MEA in the MEA Viewer and the data for the channel [24, g] in its chart.

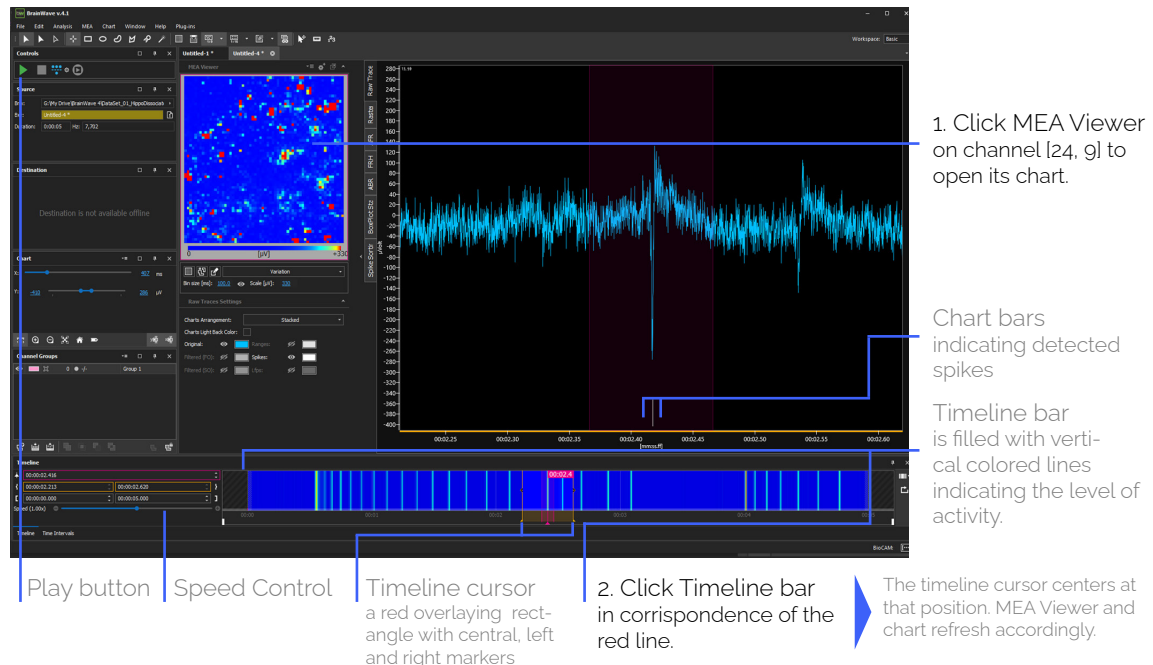

Try to move to other spatial positions in the MEA by selecting other channels in the MEA Viewer (in particular red channels) and to other time positions by clicking or dragging on the Timeline bar. To move along the time axis, you can also click Play (Stop) in order to automatically playback (stop) the data stream or alternatively you can use the Left and Right Arrow keys on the keyboard. When Play is on or you use the arrow keys, the speed at which data is refreshed is given by the Speed control under the Timeline window. Speed can also be adjusted with the Up and Down Arrow keys.

Understanding main controls affecting graphs

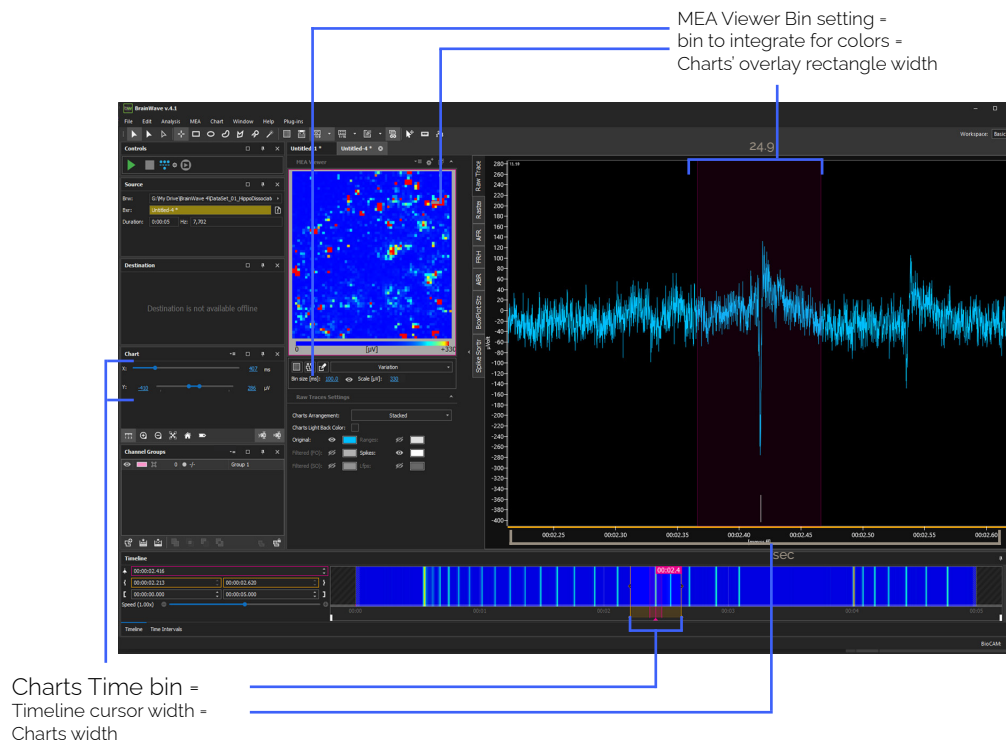


Create a group of relevant channels

You can create groups of channels according to your selection or a given statistical information. Here we create a group from a manual selection.

1. Select a group of channels of your interest by using one of the cursor tools below the main menu.
2. Click the "Create new channels group..." button at the bottom left of the Channel Groups window. A small pop-up window will appear.
3. Choose a name for the group, a labeling color, and tick the "Add Selected Channels" checkbox. Then click the "OK" button.

A group of channels will appear on the Channel Group window.

1. Cursor tools

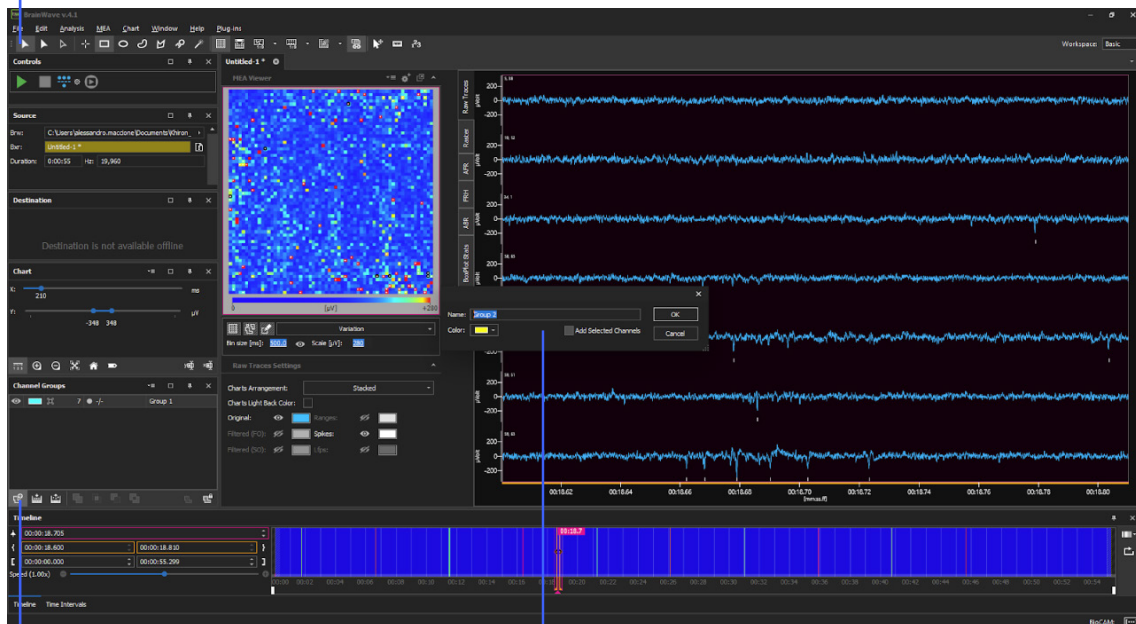

2. Create new channels group

3. Add selected channels


Visualize Instantaneous MFR

In addition to the Global MFR it might be useful to observe the instantaneous MFR, which gives the trend of the MFR computing it on successive small temporal bins. To plot the Time MFR:

1. Right-click on Channel Groups window will show the selection to select the channels belonging group. An instantaneous MFR chart is plotted by summing up together all the selected channels and binning according to the Bin size setting in the Chart window.

To zoom the MFR chart:

2. Click on the AFR window to activate the MFR chart. In this way the Charts window updates to show the controls corresponding to this chart.
3. Zoom the AFR chart either by using Chart window controls at the left or by altering the Bin value directly shown below the viewer.

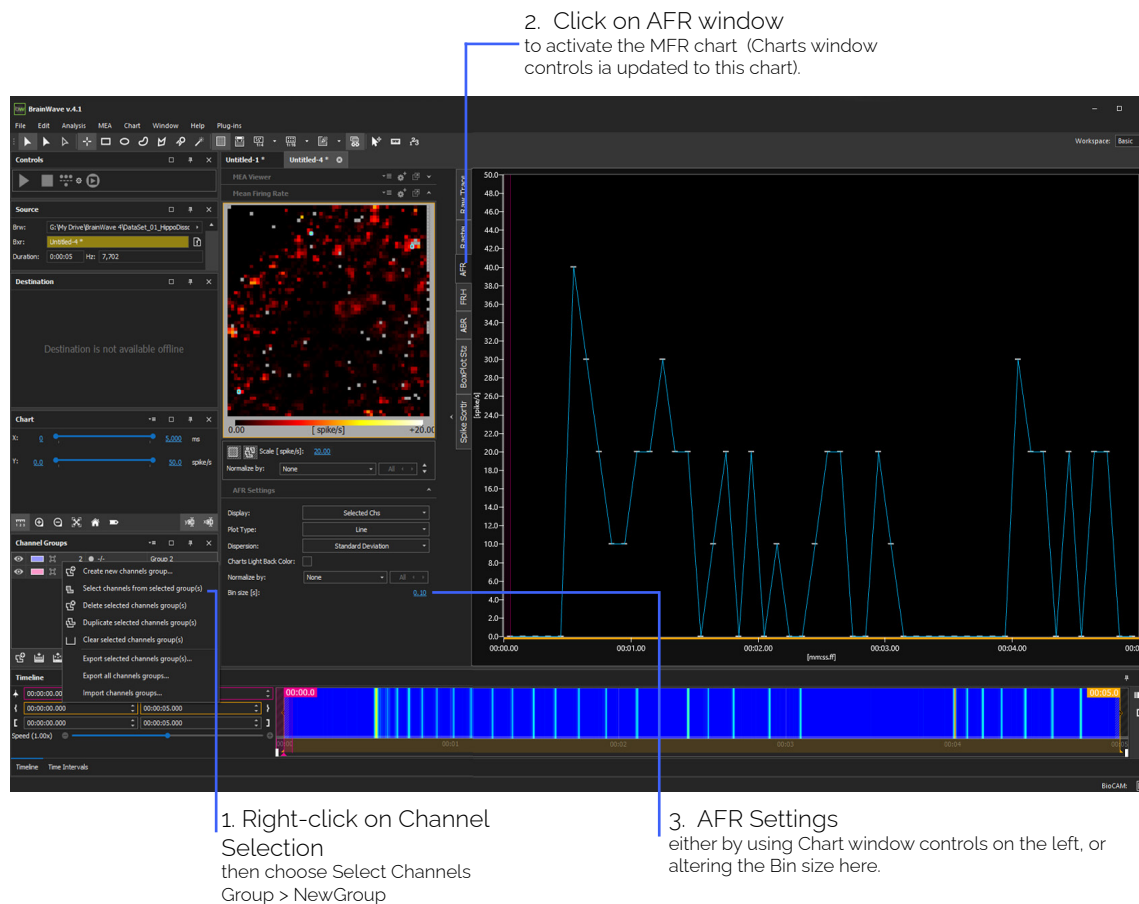


Instantaneous MFR chart is plotted by summing all the selected channels and binning according to the Bin size setting.

Visualize Raster

The time raster is another common tool used to observe activity on your data.

1. Click on Raster tab

A time raster chart is plotted.

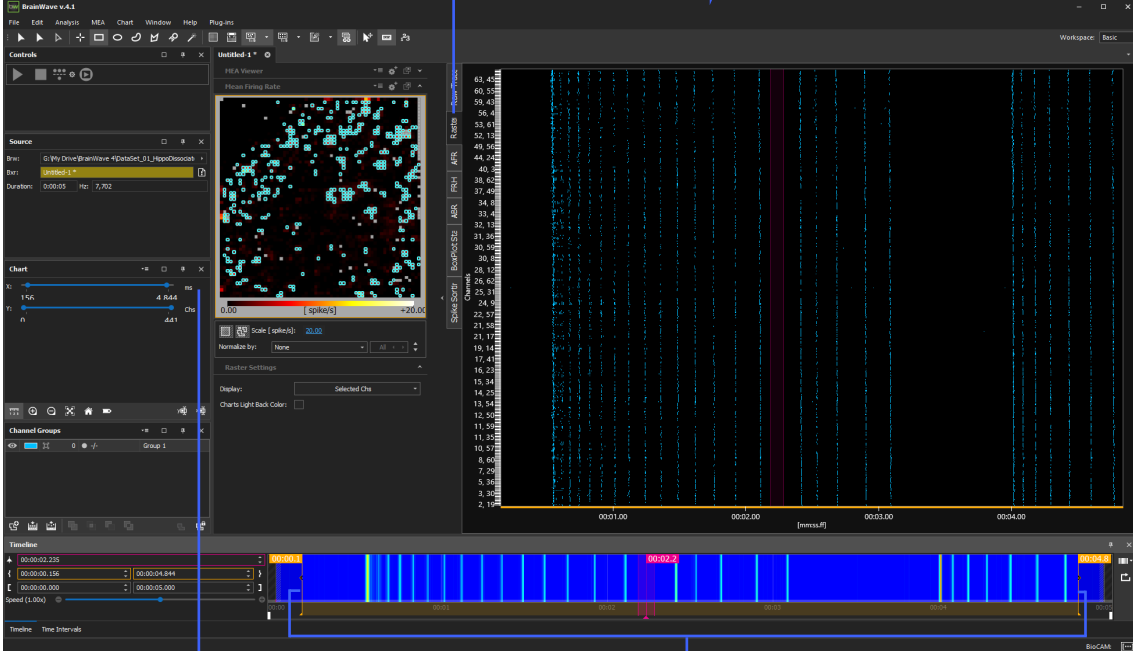

2. Decrease X axis width for the raster in order to have a more defined raster.

3. Zoom on time axis by using the Timeline bar control and moving the aside handlers.




6

Troubleshooting

Electrodes with no-signal (saturation)

Symptoms

Few electrodes in the array have either persistent or short periods with no visible signal. This effect is overall called saturation and in this section we will refer to its two possible subtypes: persistent saturation (i.e. a channel is always saturated and no signal is displayed) and jumping saturation (i.e. raw signal is affected by short periods of saturation typically of tens to hundreds of milliseconds).

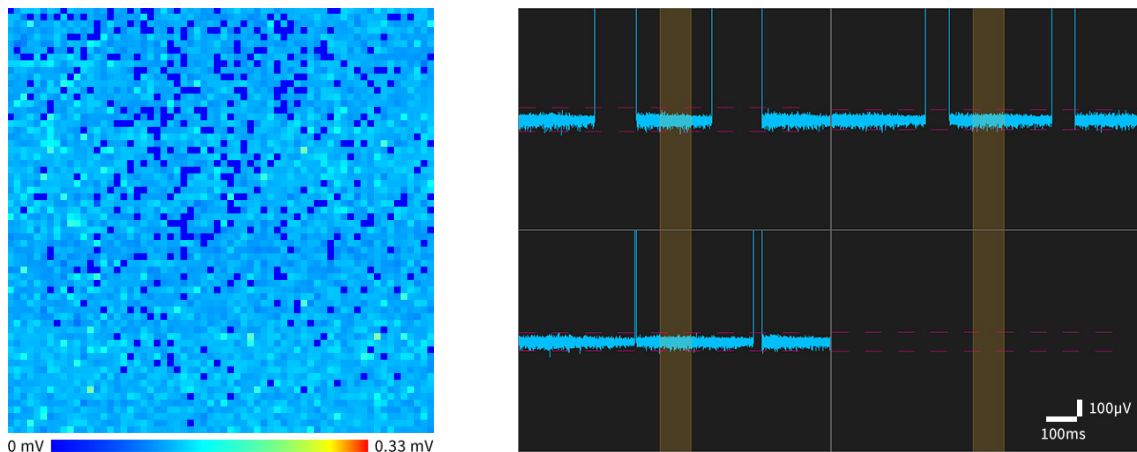

In the MEA Viewer control (left image) in BrainWave saturating electrodes show a lowest fixed color scale value (persistent saturation) or they blink toward the lowest scale value. The lowest color scale value is typically deep blue but may differ according to the chosen color scale. By displaying the raw traces of the single electrodes it is possible to see either "jumping" signals disappearing for short periods of time (right image: top-left, top-right and bottom-left traces) or no signals (right image: bottom-right trace).

Causes

Saturation arises (i) when the input signal amplitude exceeds the operational range of the amplifiers integrated inside the HD-MEA's CMOS or (ii) when the amplifiers working range is not properly set. In the first case, this could be due to invalid signal input, for instance; in the second case a wrong calibration of the HD-MEA could be the cause.

Possible scenarios and remedies


#	Scenario	Remedy
S1	Electrodes on the HD-MEA are insulated by tiny bubbles that form when adding a saline solution because of the hydrophobicity of the CMOS surface. Insulated electrodes exhibit persistent saturation because the input signal is missing. This scenario arises only when trying a new HD-MEA or a HD-MEA that has not been used for a long period. A 2x or 4x magnification lenses are enough to observe these bubbles.	Use a plastic pipette with a small tip to flow liquid on the CMOS surface in order to remove the bubbles. While doing so, be very careful not to touch the CMOS and to avoid electrodes damages. Alternatively empty the chip's reservoir, put a drop of pure etoh on the recording surface for a few seconds, rinse with water and put back in the solution

#	Scenario	Remedy
S2	CMOS is exposed to too high an environmental light source and it is not properly configured to work in such conditions. In such scenario several electrodes typically exhibit jumping saturation.	<p>Turn off the light source or cover the HD-MEA with a cap (as those that can be provided together with the BioCAM Duplex system).</p> <p>If you cannot perform this operations, or if you need light on the HD-MEA (e.g. for light stimulation purposes) configure BrainWave to make the HD-MEA working in such conditions.</p> <p>If the problem is not solved, check following scenarios and remedies.</p>
S3	This can happen for the Arena, Stimulo, and Prime chips, where the pseudo-reference electrode is dirty or it has not been cleaned for a relatively long period. Proteins and impurities can attach with time to the reference electrode and degrade the reference performances, causing few or many channels to exhibit persistent saturation.	Check section "BioCAM Duplex handling and use" P.15 how to clean the pseudo-reference electrode.
S4	HD-MEA is not properly calibrated. In such scenario some channels might show persistent saturation while others suffer from jumping saturation.	Check section "HD-MEA Calibration" P.28 to calibrate the chip.
S5	HD-MEA is either damaged or worn out.	If none of the above remedies worked you may need to change the HD-MEA.

HD-MEA not properly initialized

Symptoms

On the MEA Viewer control in BrainWave, large red horizontal stripes move up or down across the entire array.

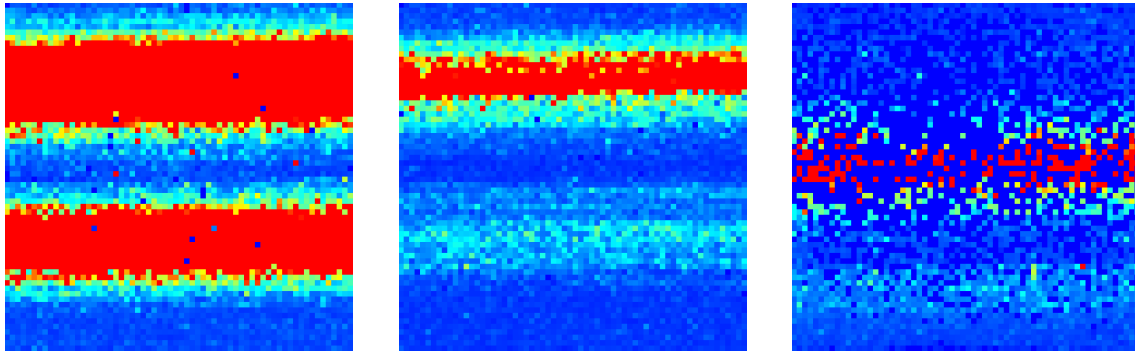

Causes

HD-MEA are monolithic CMOS-MEA chips integrating active semiconductor components (i.e. transistors) underneath and around the active area (i.e. the area exposing the sensing surface of the electrodes). Some of these active components need to be correctly biased by sending configuration settings from BrainWave to the BioCAM and the HD-MEA before starting the read out of the data. This operation is performed when starting the acquisition in BrainWave (e.g. by clicking Play button). If the connection to the HD-MEA is not in place or it is compromised, the initialization of the chip will fail.

Possible scenarios and remedies


#	Scenario	Remedy
S6	BioCAM locking system (see section "Plug the HD-MEA" P20) was in the unlocked state when acquisition was initiated.	<p>Stop the acquisition, make sure that the HD-MEA is properly inserted till the end of its bay and push the Lock button of the BioCAM Du-plex to its end position (see section "Plug the HD-MEA" P20 for more details). Make sure that the HD-MEA LED icon on the BrainWave's Right Status Bar (see section "BrainWave 4" P13) is on, signalling that the HD-MEA is properly inserted and locked. Restart the acquisition.</p> <p>If the HD-MEA LED icon on BrainWave does not turn on, see following scenarios.</p>

#	Scenario	Remedy
S7	HD-MEA contact pad is dirty or damaged.	<p>Extract the HD-MEA from the HD-MEA bay and clean all the gold contact pads on the HD-MEA by using a tissue soaked with pure ethanol (>96%).</p> <p>In case you observe some dirt on the pads you can also try to remove it by using a soft rubber. Be careful not to damage or remove the gold layer covering the electrical contacts. Then, clean again with pure ethanol and let it dry for at least one minute.</p> <p>If cleaning has no effect, try one or more different HD-MEAs to verify the problem does not lie in the BioCAM. If also other HD-MEAs exhibit the same problem, go to next scenario.</p>
S8	BioCAM DupleX locking mechanism, which is responsible to connect the HD-MEA to the BioCAM, is damaged.	<p>Double check if the locking LED shows red during the locking state. If it does, hovering over the LED should indicate the problem.</p> <p>However please contact customer support..</p>

Not working columns

Symptoms

Groups of 4 columns do not work. This typically happens on 4 columns but might also involve 8 or 12 columns and in any case a multiple of 4 columns (columns are grouped by 4 in the multiplexing logic). If you see only part of a column or a single column in a non-working conditions, then the HD-MEA is inherently damaged at that location. In most cases it is still possible to record from the remaining working electrodes.

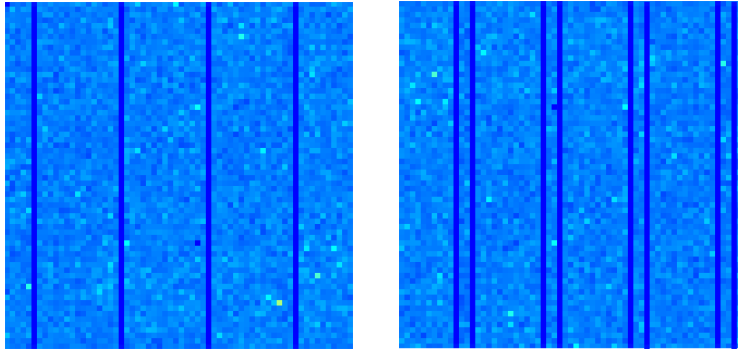

Causes

One or more of the lines that go from the HD-MEA to the BioCAM Duplex exhibit a faulty connection. This issue typically occurs between the HD-MEA contact pad(s) and the Duplex bridge module integrated inside the mechanical locking system.


Possible scenarios and remedies

Check scenarios "[S7](#)" and "[S8](#)" in previous sections.

HD-MEA with dirty surface

Symptoms

This scenario can occur when testing a HD-MEA with no biological preparation. The chip might not have been properly cleaned or has not been used for a long period. In this situation, several electrodes on the array could exhibit either higher level of noise or signals with fake peaks. In BrainWave, you typically see several spots (single electrodes) with red values on the overall array (left image). Single electrode raw traces can have either high noise levels (centre image: bottom traces are more noisy than average traces on top) or peaks at a sustained pace (right image). Most of the peaks have a shape quite different from that of a real spike, but some might still look similar.

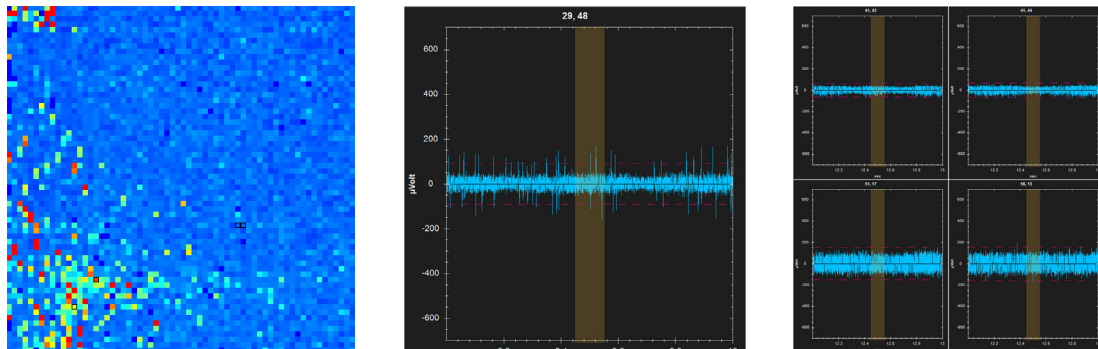

Causes

HD-MEA CMOS surface is dirty, which makes some electrodes unstable.





Possible scenarios and remedies

#	Scenario	Remedy
S9	HD-MEA electrodes are dirty of debris and/or more likely of dust particles.	<p>While the chip is connected and is filled with saline solution, start acquiring and visualizing incoming data with BrainWave to have a feed-back on whether the problem is solved.</p> <p>As a first step, try pipetting with a small plastic pipette by placing the tip roughly at 45° and very close to the CMOS central area. Be very careful not to touch the CMOS with the tip.</p> <p>If this does not help, you can also try to use a soft brush directly on the CMOS surface by gently brushing repeatedly until all the instable electrodes get recovered.</p>

Pseudo-Reference Electrode bubbles

Symptoms

This scenario occurs when bubbles have formed above the pseudo-reference electrode within the HD-MEA. This can create excess noise on the recordings.

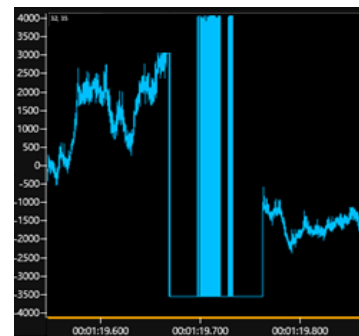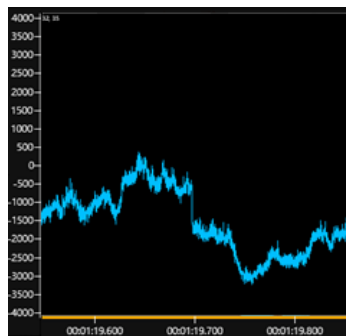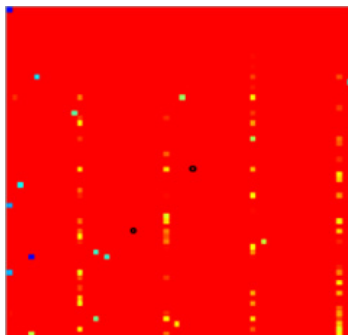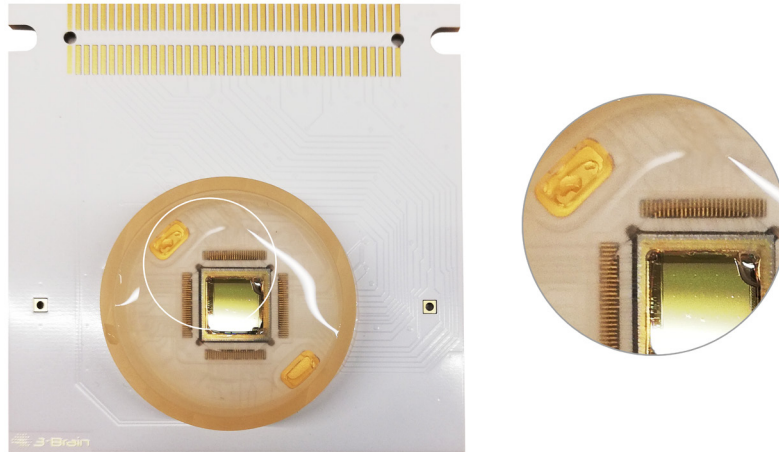

Causes

The gold pad pseudo reference is hydrophobic and some air are kept trapped at the reservoir border.

Possible scenarios and remedies

#	Scenario	Remedy
S9	Pseudo-reference electrode on HD-MEA has bubbles causing extra noise collected into the electrophysiological imaging.	Use a small plastic pipette to suck out the bubble

Contact information

If there are any further questions which you would like to personally contact us for assistance with. Please feel free to contact us through the following link:

<https://www.3brain.com/contact>

For urgent communications contact:

Phone: +41-813227086

Email: info@3brain.com

